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**Climatic, Thermostatic
and Plant Growth Chamber**

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CLIMATIC, THERMOSTATIC AND PLANT GROWTH CHAMBERS

Since 1932, Angelantoni Industrie has been designing and manufacturing climatic chambers and has become the European leader for the sector concerning environmental simulation chambers.

CLIMATIC AND THERMOSTATIC CHAMBERS FOR STABILITY TESTING

The series of EKOCH-EKOCT-EKOCTL-EKOCHL chambers introduced in this catalogue are primarily used for stability tests in the chemical-pharmaceutical sector. The small external dimensions, easy use, reduced consumption of energy and high reliability of the chambers are the main characteristics for use in research laboratories. All the models proposed comply with ICH regulations.

SOLAR LIGHT CLIMATIC AND THERMOSTATIC CHAMBERS

Solar simulation chambers enable the combination of temperature and humidity cycles with indoor lighting. They are mainly used for growing plants, insect incubation and growth, cell cultures, germination and tissues and organs.

EKOCH-EKOCT-EKOCTL-EKOCHL solar Light

Chambers provide variable lighting from 0 to 20,000 lux using cold light with hue (different hues/colours can be applied) that are fixed outside the treatment room (on both sides).

Upon request, high lighting power chambers including pre-fabricated and operator-accessible types are also available.

Angelantoni Industrie

Vista camere complete di accessori extra
View of chambers fitted with extra accessories

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STRUCTURE

Metal frame painted in grey (RAL 9006) with vertical band in blue. Back doors are provided with external lock and key. Internal double doors with toughened glass. Installed on rollers for easy handling and levelling feet included.

HEAT INSULATION

CFC-free polyurethane, (75mm average insulation thickness).

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INTERNAL CHAMBER

Stainless steel AISI 304 with internal round off corners for easy cleaning and disinfection. Height adjustable plasticized steel internal shelves.

COOLING PLANT

It is totally sealed and adopts a hermetic airtight compressor with R134a ecological refrigerant and CFC-free ODP=0. It has internal ventilation which guarantees excellent set value uniformity.

HUMIDIFICATION(EKOCH – EKOCHL models only)

Electronically-controlled low-pressure steam generator equipped with demineralised water inlet.

ADJUSTMENT AND CONTROL

Electronic digital controllers (PT 100 probe for temperature and a capacitive probe for R. H. measurements) based on the P.I.D adjustment system. Mechanical safety thermostat for internal over-heating with chamber shutdown. Alarm signal for open door, maximum temperature and internal overpressure.

Upon request, we can supply a relative temp. and humidity cycle programming system (Micro PLC control) provided with LCD screen.

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EXTRA ACCESSORIES PROVIDED ON REQUEST

- *Alarm system for : minimum and maximum temperature, min. and max. relative humidity , no mains power. The system is fitted with a buffer battery when there is no power.*
- *Electronic digital controllers for setting min. and max. temperature and relative humidity shutdown.*
- *Multipane glass heated inspection window, dimensions mm: 400 x 700 H (2 per EKOCH 1500)*
- *Additional shelves (s.s. upon request)*
- *Microprocessor graphic recorder for continuous temp. and relative humidity value recordings (including light "on-off").*
- *Single-column demineralising device for humidification system (for EKOCH – EKOCHL models only).*
- *Serial interface through RS232 or RS485 configuration for multi-drop connection to remote control system.*
- *50 mm access port hole with rubber cap to enable electrical or additional probe connections.*
- *"WinKratos" software operating on Windows 98/NT/2000/XP for controlling a chamber or for multi-drop connection.*
- *Control PC for chamber running.*
- *UV lamp with automatic shutdown switch when door is opened.*

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Angelantoni Industrie headquarters in Massa Martana (Perugia, Italy) extend over an area of 80.000 square metres (more than 16.000 square metres are covered by factories and offices).

Massa Martana is located in Umbria, a region rich in art, history and tradition.

No location could be more appropriate; **Angelantoni Industrie** S.p.A. learns from the past to better understand and anticipate the future.

This, combined with dedication and ever growing expertise, makes **Angelantoni Industrie** S.p.A. the most complete and diversified European group for advanced cold technology for industry and research. Founded in 1932, **Angelantoni Group** has today more than 600 employees working in 6 units (Italy [2], France, Germany, China, India).

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AS biomedical division also produces:

- reach-in refrigerators (+4°C and -20°C)
- plasma shock freezer
- automatic ice cubers or flakers
- -40°C and -85°C Platinum freezers
- CO₂ incubators
- chambers for pharmaceutical stability tests
- plant growth chambers (sun light)
- -152°C freezers (with compressors)
- prefabricated cold and hot chambers
- special execution refrigerators
- explosion-proof refrigerators

ISO 9001:2000

ISO 14001

AQAP 110

ANGELANTONI Industrie SpA

AS biomedical division

Loc. Cimacolle, 464 - 06056 Massa Martana (PG), Italy

Tel. (+39) 075 8955.1 - Fax (+39) 075 8955312

e-mail: biomedical@angelantoni.it

www.angelantoni.it

MORE THAN TECHNOLOGY